

Before using the accessories for the first time, read these instructions for use. Also read the operating manual of the device the accessories are used with.

## 1 Product description

### 1.1 Main illustration

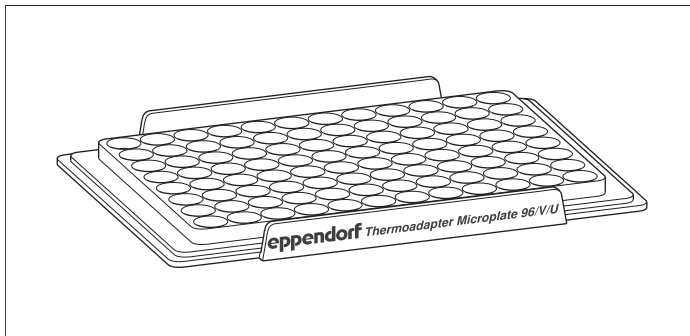


Fig. 1: Thermoadapter Microplate 96/V/U

### 1.2 Features

The Thermoadapter Microplate 96/V/U can be tempered in the *epMotion* on the Eppendorf ThermoMixer module and the thermal module. The Thermoadapter Microplate 96/V/U reduces the time in which the samples reach the target temperature.

The Thermoadapter Microplate 96/V/U can be used in any *epMotion*. The Thermoadapter Microplate 96/V/U can be used with *epBlue* from version 40.4 onwards.

The following labware can be used in the Thermoadapter Microplate 96/V/U:

- Eppendorf Microplate 96/V
- Eppendorf Microplate 96/U

The Thermoadapter Microplate 96/V/U is not transported by means of the gripper. Microplates can be placed in the adapter by means of the gripper. Microplates can be taken out of the adapter by means of the gripper.

## 2 Safety

### 2.1 Warnings for intended use



**CAUTION! Risk of burns on the hot thermoadapter.**

During operation the thermoadapter can reach temperatures of up to 96 °C.

- ▶ Do not touch the hot thermoadapter.
- ▶ Wait until the thermoadapter has completely cooled down.

## 3 Ordering information

Order no. (International)	Description
5075 751.577	<b>Thermoadapter Microplate 96/V/U</b> for Microplate 96/V/U
0030 601.300	<b>Microplate 96/V-PP</b> 80 plates, wells clear, white border color PCR clean
0030 602.307	Sterile
0030 603.303	DNA LoBind
0030 601.670	<b>Microplate 96/V-PP</b> 80 plates, wells white, gray border color PCR clean
0030 601.904	<b>Microplate 96/V-PP</b> 80 plates, wells black, white border color PCR clean
0030 601.203	<b>Microplate 96/U-PP</b> 80 plates, wells clear, white border color PCR clean
0030 602.200	Sterile
0030 601.572	<b>Microplate 96/U-PP</b> 80 plates, wells white, gray border color PCR clean
0030 601.807	<b>Microplate 96/U-PP</b> 80 plates, wells black, white border color PCR clean

Your local distributor: [www.eppendorf.com/contact](http://www.eppendorf.com/contact)

Eppendorf AG · 22331 Hamburg · Germany  
[eppendorf@eppendorf.com](mailto:eppendorf@eppendorf.com) · [www.eppendorf.com](http://www.eppendorf.com)